# COMET4

[www.test4all.es](http://www.test4all.es)

**ZEULAB**  
Making Testing Easier

ZEULAB, S.L.  
C/ Bari, 25 dpdo. • 50197 Zaragoza (SPAIN)  
Tel.: +34 976 731 533  
[info@zeulab.com](mailto:info@zeulab.com) • [www.zeulab.com](http://www.zeulab.com)

Comet4 is an automatic device for the detection of antibiotics, controlled from the mobile phone or tablet. Results are displayed in real time wherever you are and are stored in the Test4all cloud for consultation whenever you need them. Comet4 required the Eclipse Farm 4G kit for analysis.

## **COMET4 CONNECTIVITY**

The smartphone communicates with Comet4 via bluetooth to launch the assays. If Comet4 has no Wi-Fi communication, results will be uploaded to the cloud when the device and smartphone connect. Comet4 sends the data to the Test4all cloud via Wi-Fi and provides results information in real time wherever you are and without the need to be near the device.

**Offline Mode.** If there is no internet coverage, the device can be used in offline mode. The results can be seen in the app when the smartphone and Comet4 are connected by bluetooth. When Comet4 has access to the internet using Wi-Fi or mobile coverage, the results will be synchronized in the cloud and emails will be sent.

## **ANDROID / IOS MODE**

The user can switch between Android mode and iOS mode, depending on the Smartphone operating system to be used with Comet4.

When connecting Comet4, either with cable or battery: the light flashes three times in green if Comet4 is in Android mode or three times in white if it is in iOS mode.

To change the mode:

1. Hold the button down for 10s, visually there are 10 flashes of light in red or green depending on whether it was heating or ready for use. It is not possible to change the mode if a test is in progress (blue light).
2. If the process has been completed successfully Comet4 will restart and show the new mode with the light code described above.

## INDICATIONS FOR USE AND BATTERY

Comet4 can be used connected to the mains or to the battery module. The device only has a touch button that responds quickly to soft touches.

Above the touch power button, a light indicates the status of the device.

Fixed lights:

- Red light: the device is warming up to reach the analysis temperature. It takes approximately 10 minutes.
- Green light: The device is ready to start a scan.
- Blue light: test in progress. Comet4 cannot be turned off until the test ends or is canceled.
- Pink light: Comet4 battery is charging.

Flashing lights:

- 3 green flashes when turning on the equipment. Comet4 is configured for use in Android mode.
- 3 white flashes when turning on the equipment. Comet4 is configured for use in iOS mode.
- Yellow flashes: Comet4 is updating the device firmware.

### Auto power off

For security, Comet4 will shut down if not used after 1 hour.

### Battery charge

The battery is charged when it is coupled to the Comet4 and it is turned off and connected to the mains. A pink light indicates that the battery is charging. After a full charge, approximately 12 hours, the light turns off. If Comet4 is on (red or green light), the battery does not charge.

In a continuous use of the device, it is recommended to have a full charge to launch the last test of the day with the equipment plugged in and with the battery coupled. One hour after the end of the test, the equipment will automatically turn off and the battery will begin charging.

## **Battery life**

The battery run-time is approximately 8 hours.

## **Comet4 reset**

To reset the device, disconnect the power (plug or battery), connect the device to the mains and turn on the equipment with the touch button.

## **COMET4 APP**

The Comet4 application (app) is for a comfortable handling of the device from your smartphone or tablet.

The Comet4 app can be downloaded from the Google Play Store (Android) and from the Apple App Store (iOS). Enter "Comet4 Zeulab" for searching the app.

## **TEST4ALL ACCOUNT**

The Test4all account allows the management of devices, users and results. There are two categories of users: "Admin" and "User":

- "Admin" allows registering and modifying devices and users and accessing all the registered results and devices of the company.
- "User" can only use the assigned devices and see the results to which authorization has been given.

In the "Results" section, you can see the results of each of the registered teams and download all the data in CSV format.

### **Create a Test4all account**

Account creation can be accessed from [www.test4all.es](http://www.test4all.es) by clicking on "Sign up" or from the app by selecting "Create account".

Several fields must be completed and the device key, that appears on your Comet4 certificate, must be entered. Click on "Confirm account" to complete the process.

## USE OF COMET4

In order to use Comet4, it requires to connect the device and the smartphone or tablet.

1. Turn on Comet4, open the app and log in with your Test4all account.
2. Tap "Connect" and select your Comet4 device from the list. Wait for it to connect.

### Run a test

Comet4 must be at the test temperature (63°C). It takes approximately 10 minutes for the device to reach that temperature and during this process it will show a red light. When it has reached it, it will show a green light and the device is ready to start the test.

1. Tap on "Run test".
2. Select the position (1 to 4) where the sample will be placed.
3. Tap on "Scan new batch code" and scan the QR provided in the Eclipse certificate when a new batch is started. If you are using the same batch from previous assays, tap on "Use same batch code".
4. Tap on "Test Details" to enter the identification name of the sample and tap on "Save" (this step is optional).
5. Cut an Eclipse tube with scissors (take care not to damage the foil on the remaining tubes). Add 100 µL of milk with the provided pipette, seal with the film and place it into the chosen position of Comet4 (1-4).
6. Select "Run test".
7. Comet4 will stop automatically and report the results.

At Test4all.es you can monitor the progress of the test at any time. If Comet4 and the smartphone are connected by bluetooth, the app will also show this progress in the "Status" section.

## RESULTS

Once the test is finished, if Comet4 has Wi-Fi, the results are uploaded to the Test4all.es cloud where they can be consulted at any time, and emails are automatically sent to the configured accounts.

From Test4all.es you can download the results to a ".csv" file to open them with Excel.

When Comet4 and the smartphone are connected by bluetooth, the app will show the last 100 results carried out with that device on the results screen. Clicking on a result you can see the complete information.

This information can be shared to another email from the "Share" button.

If the test has been carried out in "Offline" mode, results can be seen in the app when the smartphone and Comet4 are connected by bluetooth. When Comet4 has internet access using your wi-fi or smartphone internet coverage, the results will be synchronized in the cloud and emails will be sent.

## SETTINGS

In this section of the app, you can do different actions. The app must be connected to Comet4.

**Language:** allows you to select the language.

**User:** allows you to close the user session. The next time the app opens, it will ask for the username and password.

**WiFi:** allows you to configure the WiFi of the Comet4 equipment.

1. Select "Scan" to find a Wi-Fi network. Depending on the connection it may be necessary to touch "Scan" several times
2. Select the desired Wi-Fi network, enter the password and click "Test" to check the connection. When it says "Success", tap "Save" to save.

**About:** View the firmware version loaded on the device and update it. Tap "Update" and confirm that you want to update the device in the question "Do you want to update the firmware of the device?". If there is a higher

version, Comet4 will be updated. During the update process the light blinks yellow. After the update is complete, Comet4 restarts. Entering this section again you can see the new firmware version installed.

In this section you can also see the version of the Comet4 application.

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